## EV-Perm™ Permeabilization Pretreatment Kit for Exosome Membrane

Please visit our website through URL below and enter product code (294-85701) to learn more about product features and instructions.

<https://labchem-wako.fujifilm.com>

### 【Product Information】

Code No.	Product Name	Volume	Storage
294-85701	EV-Perm™ Permeabilization Pretreatment Kit for Exosome Membrane	1 kit	Frozen (-20℃)

### 【Kit Contents】

Reagent	Volume	Storage
Reagent A (20×)	1×500 μL	Frozen (-20℃)
Reagent B (10×)	1×5 mL	Frozen (-20℃) before thawing Refrigeration (2–10℃) after thawing
Reagent C (100×)	1×500 μL	Frozen (-20℃) before thawing Refrigeration (2–10℃) after thawing

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EV-Perm™ Permeabilization Pretreatment Kit for Exosome Membrane (EV-Perm™) enables the detection of internal extracellular vesicle proteins when PS Capture™ Exosome ELISA Kit series or PS Capture™ Exosome Flow Cytometry Kit are used in combination with EV-Perm™.

This product is research use only.

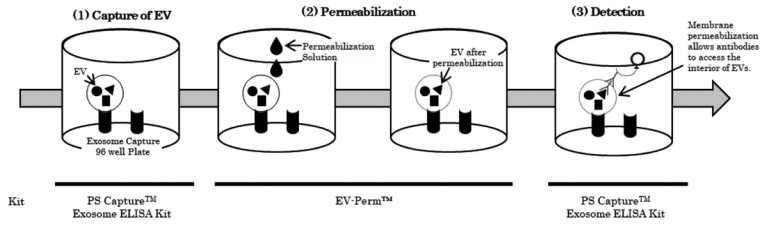
### 【Attention】

When using EV-Perm™, it is necessary to combine it with one of the following.


Code No.	Product Name
297-79201	PS Capture™ Exosome ELISA Kit (Anti Mouse IgG POD)
298-80601	PS Capture™ Exosome ELISA Kit (Streptavidin HRP)
297-79701	PS Capture™ Exosome Flow Cytometry Kit

**[Procedure for detection using EV-Perm™]**

■ If you use EV-Perm™ in conjunction with the PS Capture™ Exosome ELISA Kit series

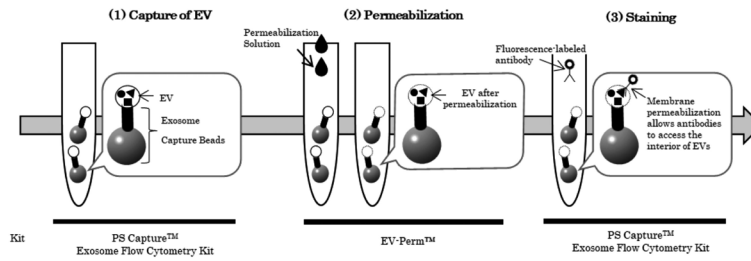■ If you use EV-Perm™ in conjunction with the PS Capture™ Flow Cytometry Kit

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**FUJIFILM Wako Chemicals U.S.A. Corporation**
U.S.A.
**FUJIFILM Wako Chemicals Europe GmbH**