

<How to replace the Pressure Adapter>




Plug

- (1) Turn off the power.

Disconnect the plug of the power cable from the outlet.

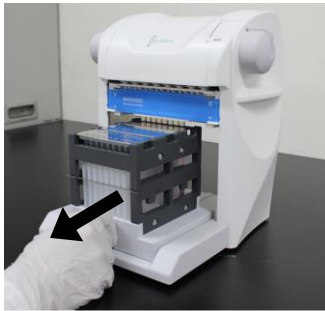

- (2) Pull out the Cartridge Holder and Collection/Waste Tube Holder.

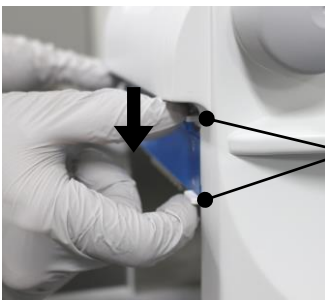

Knobs

- (3) Hold the pressing type knobs of both sides of the Mini480 Pressure Adapter as shown in the figure, and then press the upper knobs in the direction of the arrow.



-
- (4) While pressing the upper knobs, take the upper protruding parts out of the QuickGene-Mini480 as shown.
-

Note:

Do not use excessive force otherwise it may cause the Pressure Adapter failures.



Upper protruding parts

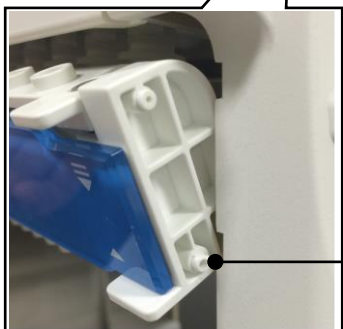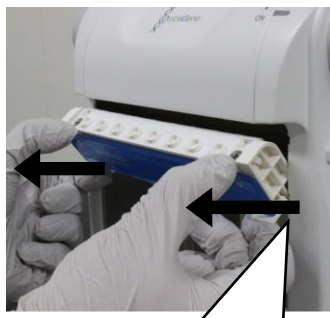


Lower protruding parts


- (5) Take out the lower protruding parts of Pressure Adapter in the direction as shown.



- (6) Hold the pressing type knobs of both sides of the Mini8L Pressure Adapter and put in the lower protruding parts of Pressure Adapter in the direction as shown.

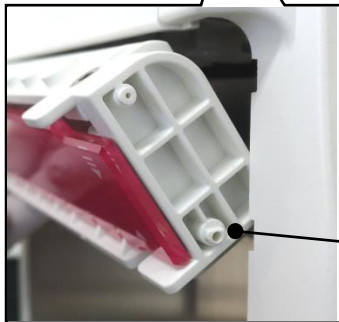


Lower protruding parts

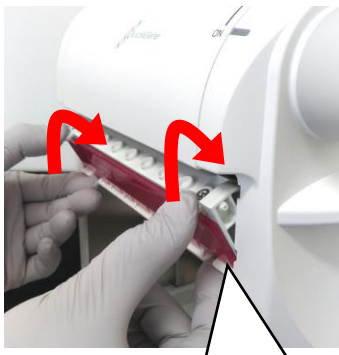


- (7) While pressing the both side of knobs, put the upper protruding parts in of the QuickGene-Mini480 as shown.

Note:

Do not use excessive force otherwise it may cause the Pressure Adapter failures.

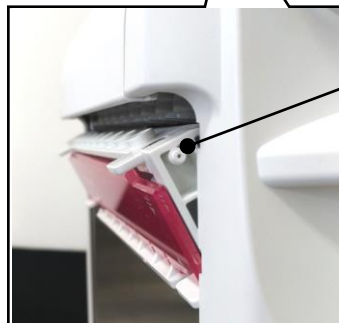

Upper protruding parts

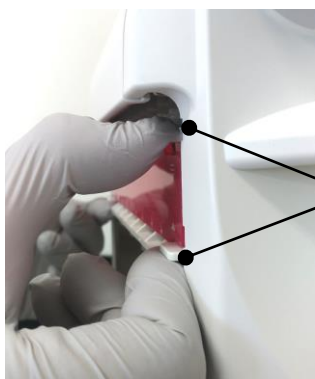

Knobs

- (8) Set the Mini8L Pressure Adapter as shown in the figure, and then release knobs.