

FOR RESEARCH USE ONLY www.genovis.com STORE AT



# **Smart**Enzymes™





### INSTRUCTIONS FOR PRODUCT

FucosEXO<sup>™</sup> 2000 units (G1-FM1-020) Digestion of up to 2 mg glycoprotein

Last revised Mar 2021

## QUICK GUIDE



 Reconstitute FucosEX0 in 50 µl ddH<sub>2</sub>0 to a concentration of 40 units/µl



Add 1 unit FucosEX0 / 1 µg glycoprotein



• Incubate for 1 to 2 h at 37°C



## PRODUCT DESCRIPTION

FucosEXO is a mix of a-fucosidases for efficient removal of a1-2, a1-3 and a1-4-linked fucose residues on *N*- and *O*-glycosylated proteins or oligosaccharides. FucosEXO hydrolyzes glycoproteins under native conditions and displays a high activity in a pH range from 6 to 8. The enzymes in FucosEXO are expressed in *E. coli*. FucosEXO is composed of two fucosidases modified with His-tags and with the molecular weights of 87 kDa and 64 kDa.

### **Unit Definition**

One unit of FucosEXO hydrolyzes fucoses from ≥90% of 11 nmol 2'-fucosyllactose and 4.5 nmol 3-fucosyllactose when incubated in 20 mM Tris pH 6.8 at 37°C for 30 min.



### **Content and Storage**

FucosEXO is supplied lyophilized in TBS pH 7.6, with no preservatives added.

FucosEXO is shipped cold and should be stored at -20°C upon arrival.

After reconstitution FucosEXO is stable for at least 1 month at +4-8°C.

FucosEXO is for R&D use only.

### DETAILED PROTOCOL

#### Additional Materials Required

• Reaction buffer1: 20 mM Tris pH 6.8

#### Sample Preparation

 Prepare the glycoprotein of interest in the reaction buffer at a concentration of 0.5-5.0 mg/ml.

#### **1** Prepare FucosEXO<sup>™</sup>

 Reconstitute FucosEXO in 50µl ddH<sub>2</sub>O to a concentration of 40 units/µl.

#### 2 Add FucosEXO<sup>™</sup>

Add 1 unit FucosEXO / 1 µg glycoprotein<sup>2</sup>.

#### **3** Digestion

Incubate for 1 h to 2 h at 37°C

Optimization of enzyme concentrations and incubation time may be needed depending on the substrate.

#### Notes

- FucosEXO displays high activity in buffers with pH values from 6 to 8 and over a wide range of ionic strength (0-500 mM NaCl). Some optimizations might be required if a buffer other than the recommended reaction buffer is used.
- A higher enzyme concentration may increase digestion efficiency of individual glycoproteins. This requires optimization.



### **Quality Control**

FucosEXO is tested to meet the specifications and lot-to-lot consistency.

FucosEXO is tested for the absence of microbial contamination using blood agar plates, Sabouraud dextrose agar plates and fluid thioglycollate medium.

### **Related Products**

SialEXO<sup>®</sup> For complete removal of  $\alpha$ 2-3,  $\alpha$ 2-6 and  $\alpha$ 2-8 linked sialic acids.

#### GalactEXO<sup>™</sup>

For complete removal of  $\beta$ 1-3 and  $\beta$ 1-4 linked galactoses.

#### Legal and Disclaimers

All rights reserved. Genovis products may be covered by one or more patents, trademarks and copyrights owned or controlled by Genovis AB.

For more information about commercial rights, please contact Genovis team at licensing@genovis.com

Genovis products are intended for research use only. They are not intended to be used for therapeutic or diagnostic purposes in humans or animals.

© Genovis AB



#### **USA & Canada**

Genovis Inc. 245 First Street, Suite 1800 Cambridge, MA 02142 USA

Customer service: 1-617-444-8421 Order phone (toll free): 1-855-782-0084 Order fax: 1-858-524-3006 Email: orders.us@genovis.com

#### **EMEA & Asia**

Genovis AB Box 790 SE-220 07 Lund Sweden

Customer service: +4646101230 Order phone: +4646101230 Order fax: +4646128020 Email: order@genovis.com