




GSEV&ASEV meeting

Characteristics on PS Affinity for Isolation and Detection of EVs: Advantages Clarified from Comparison with Conventional Methods

Key word: extracellular vesicles, exosome, microvesicle, phosphatidylserine

Sadamu Ozaki 尾﨑 真夢 Strategy Planning & Coordination Department Laboratory and Specialty Chemicals Division FUJIFILM Wako Pure Chemical Corporation

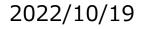

Today's Outline

- **1.** Introduction on PS Affinity
- 2. Comparing PS Affinity with Ultracentrifugation and SEC
- 3. Reference Data -Proteomic Analysis -Density Distribution
- 4. Conclusion

1. Introduction on PS Affinity

Introduction on PS Affinity Method FUJiFILM







MagCapture[™] Exosome Isolation Kit PS Launched on Dec in 2015


- Collaborated with Prof. Rikinari Hanayama. (WPI Nano Life Science Institute, Kanazawa Uni.)
- Developed in 2015
- Isolating and Detecting EVs (e.g. Exosome etc.)
- Completely different from conventional methods
- Introducing highly-pure extracellular vesicles compared with conventional methods



OPEN A novel affinity-based method for the isolation of highly purified extracellular vesicles

Nakai et al., Sci Rep. 2016 Sep 23;6

EV Research Products using PS Affinity

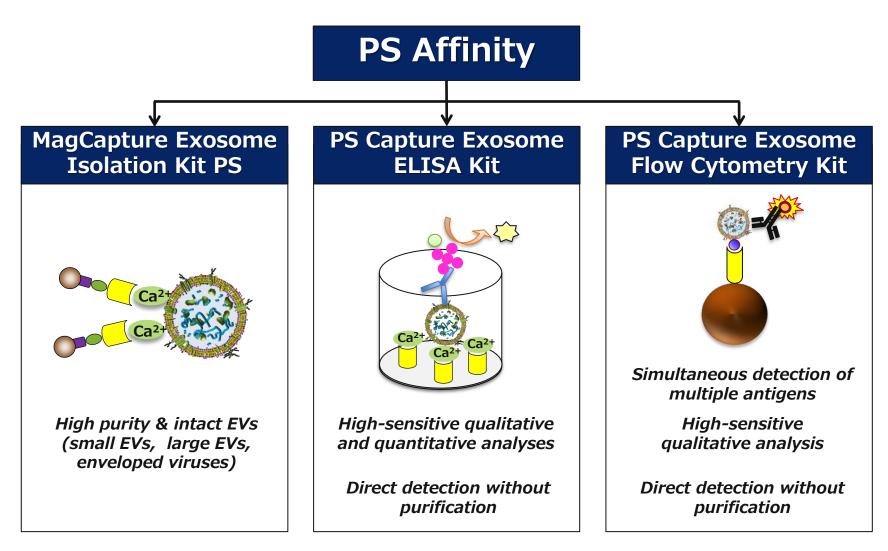


Easy operation and high reproducibility

Characteristics of PS Affinity Method

FUJIFILM

Scheme of phosphatidylserine (PS) affinity-based method

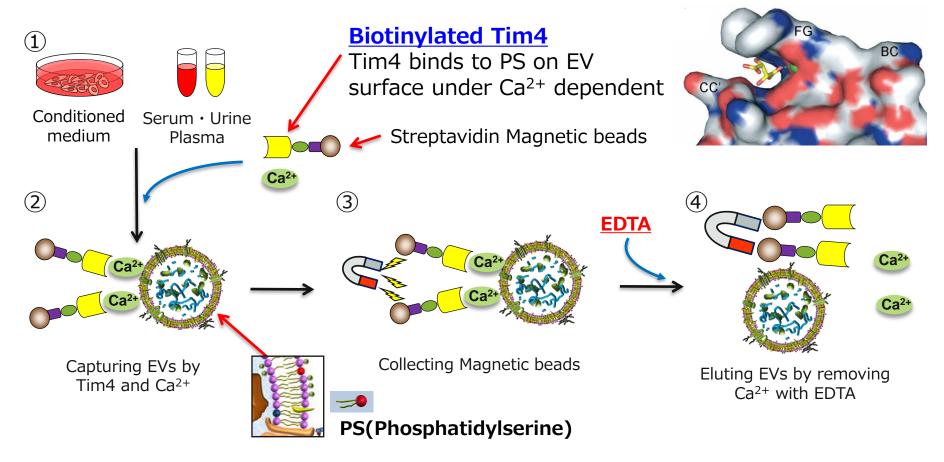




Wako

<u>A novel affinity purification method is capable of purification of highly pure</u> and intact extracellular vesicles

Investigating Superiority of PS Affinity

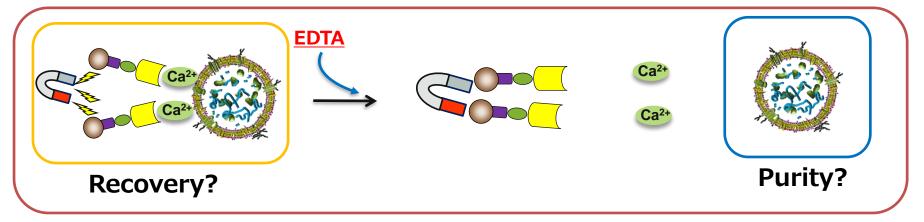

Validity?

Evaluation Points of PS affinity method

- 1. Recovery amount
- 2. Purity
- 3. Validity as method

For proving superiority in terms of the above, we carried out comparison between conventional methods and commercial kit.

2. Comparing PS Affinity with Ultracentrifugation


- 1. Appearance
- 2. Purity
- 3. Recovery Amount and Recovery Efficiency

Comparing PS Affinity with UC


- in terms of appearance by TEM analysis -

■ sEVs in 10K sup of COLO201 cells were isolated by each method, followed by transmission electron microscope (TEM)

PS Affinity

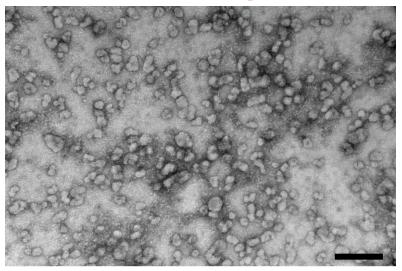

Ultracentrifugation

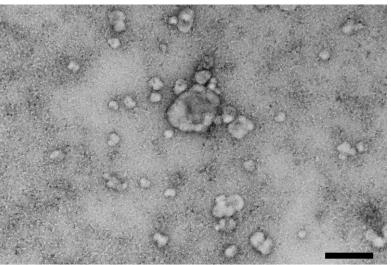

x 100,000 Bar : 200nm


FUJ!FILM



PS Affinity isolated many particles of around 50-100 nm in size. Numerous small EVs can be seen.

Comparing PS Affinity with UC

- in terms of purity of small EVs by biochemical analysis -

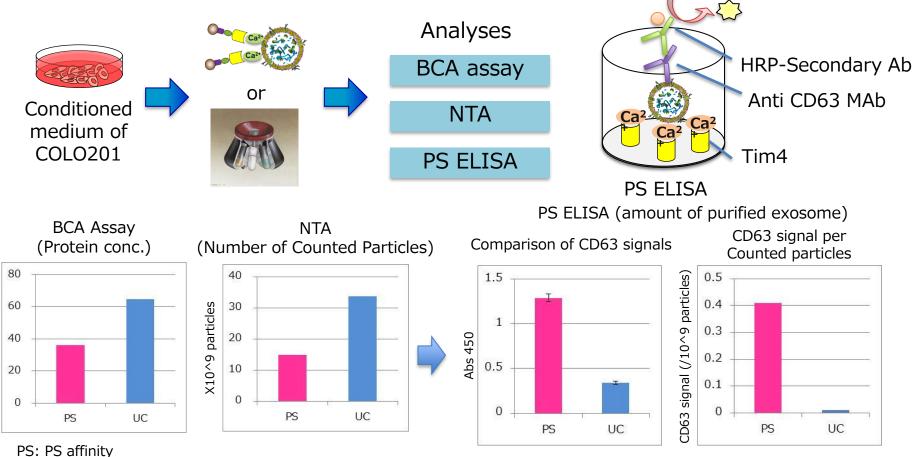


UC: Ultracentrifugation

Wako

²rotein Concentration (ng/µl)

PS affinity method enables the isolation of highly pure small EVs.

Comparing PS Affinity with UC

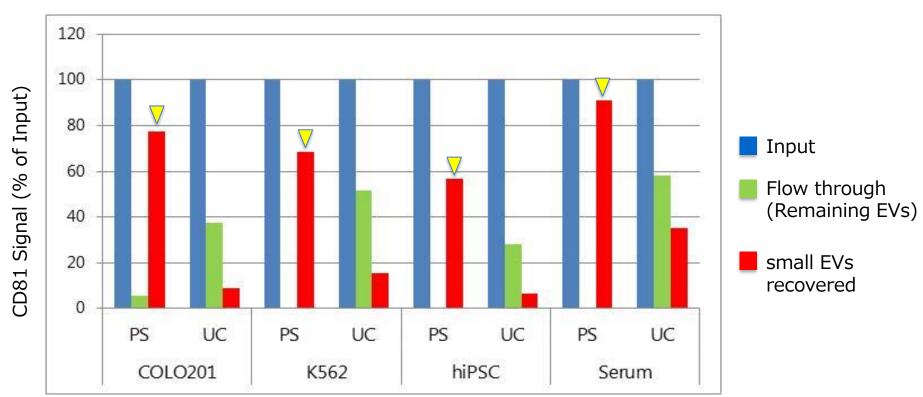

- in terms of recovery efficiency by PS ELISA -



These results indicates that PS affinity can recover sEVs derived from various cell lines and serum more efficiently than ultracentrifugation.

3. Reference Data

- **1. Proteomic Analysis**
- 2. Density Distribution

Proteomic Analysis

EV markers

Wako

12


Top 15 proteins identified in the isolated sEVs

	Tim4 method	Ultracentrifugation	PEG precipitation
1	Heat shock cognate 71 kDa protein	DNA-PK catalytic subunit	Complement C3
2	Annexin A6	Transferrin receptor protein 1	Alpha-2-macroglobulin
3	Transferrin receptor protein 1	Serum albumin	Fibronectin
4	V-type proton ATPase subunit A	ATP-dependent RNA helicase A	Serum albumin
5	Flotillin-2	Tubulin beta-5 chain	Thrombospondin-1
6	Programmed cell death 6	Heat shock cognate 71 kDa protein	Complement C4
7	4F2 cell-surface antigen heavy chain	Fatty acid synthase	Alpha-1-antiproteinase
8	Annexin A1	4F2 cell-surface antigen heavy chain	Apolipoprotein B-100
9	Kinase D-interacting substrate	U5 small nuclear RNP helicase	Hemoglobin fetal subunit beta
10	Annexin A2	Tubulin beta-4B chain	Tubulin beta-5 chain
11	Flotillin-1	Ribonucleoprotein M	Fatty acid synthase
12	V-type proton ATPase subunit B	Hemoglobin fetal subunit beta	Adiponectin
13	Annexin A11	Clathrin heavy chain 1	Fibulin-1
14	Annexin A7	Fibronectin	Complement C4A
15	Syntenin-1	Tubulin alpha-1B chain	Complement C7



Gray columns: bovine proteins from FBS

Density distribution of isolated sEVs FUJIFILM

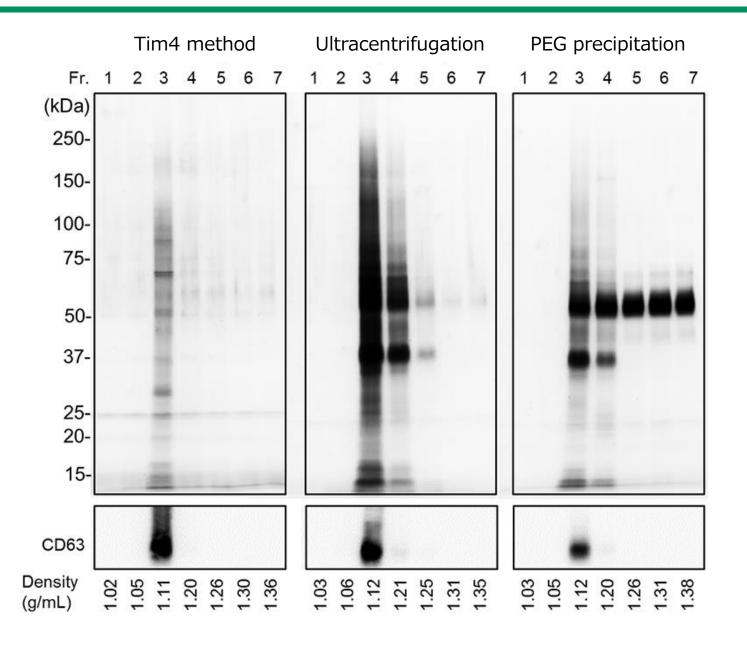


Wako

4. Conclusion

Conclusion

FUJIFILM

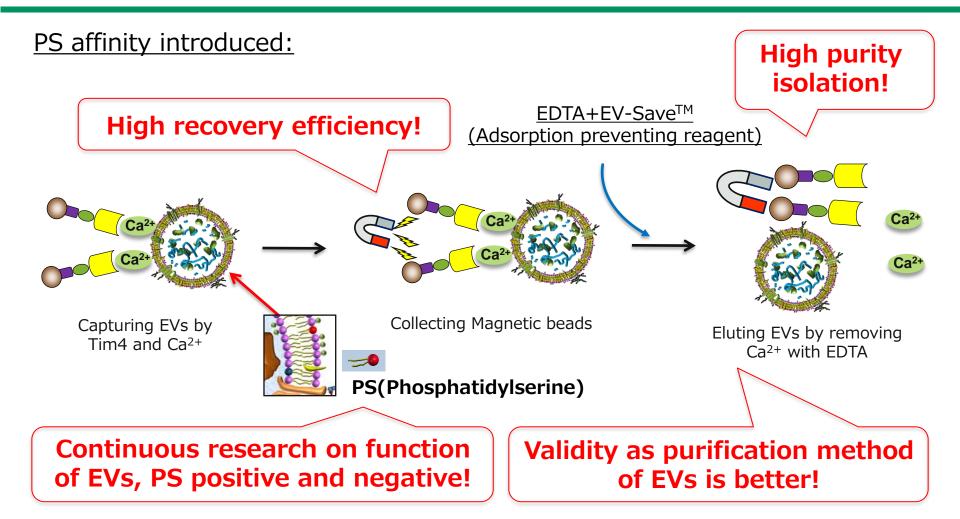

PS Affinity exhibited the superiority to conventional method such as ultracentrifugation and also presented new challenge!

~Additional Information~ Do you know the loss and adsorption of EVs to labwares?


Features

Strongly suppressing adsorption of EVs to laboratory tools

EV-Save[™] Extracellular Vesicle Blocking Reagent


- Simple operation just to add to the sample
- Cryoprotective effect to EVs

Tube Transfer

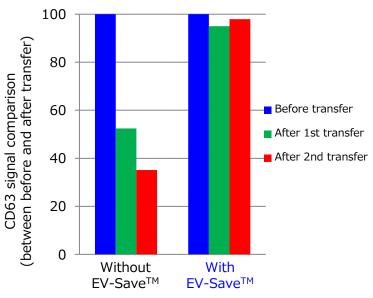

Ultrafiltration


Description	Package Size	Catalog No.	Storage	
EV-Save [™] Extracellular Vesicle Blocking Reagent	1 mL	058-09261	Keep at -20℃.	17

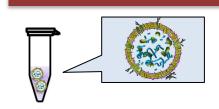

Save EVs

Cryoprotective effect by EV-Save™

Repeating freeze & thaw

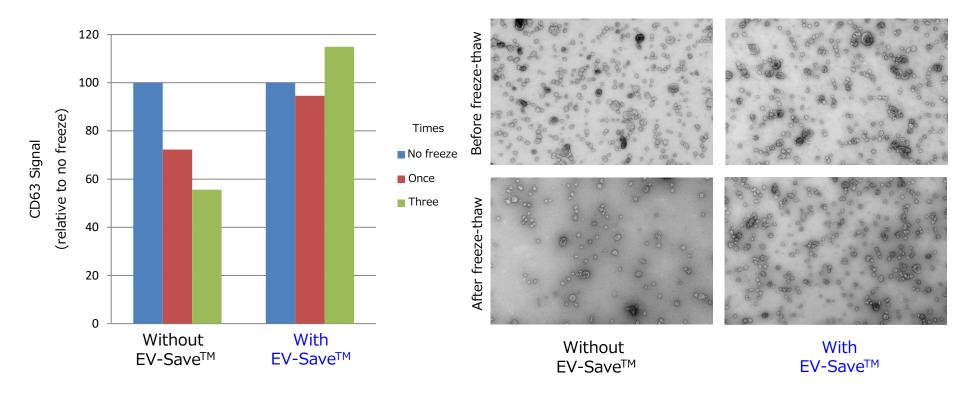

[Result]

Although freeze-thaw reduced CD63 signal, such reduction was suppressed by EV-Save[™] (A). Furthermore, TEM results (B) indicated that freezing and thawing caused a marked decrease in the number of particles, but the EV-Save[™] suppress it.

FUJ:FILM Value from Innovation