

(90 × 190mm Size)

FUJIFILM

Wako

Code No. 043-31381 (1 g)

049-31383 (5 g)

Deoxyribonucleic Acid, from Salmon Sperm

デオキシリボ核酸, サケ精液由来

This product is the sheared DNA from salmon sperm. It is used to prevent non-specific binding of the probe to the membrane in Southern and Northern Hybridization. This product should be sonicated, extracted with phenol/chloroform and precipitated with ethanol before using.

[**Appearance**] Lyophilized

[**Size of DNA**] Approximately 500bp~1,000bp

[**Storage**] 2~10℃ · 遮光保存

[**Package**]

Code No.	Package
043-31381	1 g
049-31383	5 g

FUJIFILM Wako Pure Chemical Corporation

1-2, Doshomachi 3-Chome, Chuo-Ku, Osaka 540-8605, Japan
Telephone : + 81-6-6203-3741
Facsimile : + 81-6-6201-5964
<http://www.wako-chem.co.jp>

FUJIFILM Wako Chemicals U.S.A. Corporation

1600 Bellwood Road
Richmond, VA 23237
U.S.A.
Telephone : + 1-804-271-7677
Facsimile : + 1-804-271-7791
<http://www.wakousa.com>

FUJIFILM Wako Chemicals Europe GmbH

Fuggerstrasse 12
D-41468 Neuss
Germany
Telephone : + 49-2131-311-0
Facsimile : + 49-2131-311100
<http://www.wako-chemicals.de>

1801KA1