(90 × 210mm Size)

FUJIFILM

Wako

Code No. 193-18455

For Tissue Optical Clearing Reagent
SCALEVIEW-A2

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Dr. Atsushi Miyawaki et al. developed a water-based optical clearing reagent, Scale, which clears fixed biological samples without quenching fluorescent protein.

SCALEVIEW-A2, bodily tissues can be cleared of their opacity to provide a clearly defined image.

【Additional required materials】

Reagents :

- 1) Paraformaldehyde (Code No. 160-16061)
- 2) 1 × PBS(-) (Code No. 164-25511)
- 3) Sucrose (Code No. 194-00011)
- 4) Immersion solution

Please use the immersion solution recommended by microscope manufactures.

Equipments :

- 1) 50 mL conical centrifuge tube
- 2) Seesaw shaker (for fragile samples)
- 3) Recommended fluorescence microscopes and objectives for SCALEVIEW-A2 treated samples

1. Multiphoton microscope : OLYMPUS, model : FVMPE-RS

● Multiphoton dedicated objectives :

- OLYMPUS, model : XLPLN25XSVMP : 25 ×, NA = 1.0, WD = 4 mm
- OLYMPUS, model : XLSLPLN25XSVMP : 25 ×, NA = 0.95, WD = 8 mm
- OLYMPUS, model : XLPLN10XSVMP : 10 ×, NA = 0.6, WD = 8 mm

● Objectives :

- OLYMPUS, model : UPLSAPO30XSIR : 30 ×, NA = 1.05, WD = 0.8 mm

2. Confocal microscope : OLYMPUS, model : FV3000

● Objectives :

- OLYMPUS, model : UPLSAPO30XS : 30 ×, NA = 1.05, WD = 0.8 mm
- OLYMPUS, model : UPLSAPO40XS : 40 ×, NA = 1.25, WD = 0.3 mm
- OLYMPUS, model : UPLSAPO60XS : 60 ×, NA = 1.30, WD = 0.3 mm
- OLYMPUS, model : UPLSAPO100XS : 100 ×, NA = 1.35, WD = 0.2 mm

【Procedure】

A case for making mouse brain transparent according to reference (1).

- 1) Transcardially perfuse an anesthetised mouse with 4 % paraformaldehyde (PFA)/PBS (pH 7.5~8.0).
- 2) Remove the whole brain and subject it to post-fixation in 4 % PFA/PBS at 4 °C for 10 hrs and cryo-protection in 20% sucrose/PBS at 4 °C for 24 hrs.
- 3) Embed the sample in OCT compound and freeze it with liquid nitrogen.
- 4) Thaw and rinse the sample in PBS, and fix it again with 4 % PFA/PBS for 20 min at room temperature.
- 5) Clear the sample by incubation in SCALEVIEW-A2 at room temperature.

More than 30 mL of SCALEVIEW-A2 is required for an adult mouse brain.

A full week of incubation may be necessary for transparency. During incubation, stir slowly using an orbital shaker.

Exchanging SCALEVIEW-A2 each day accelerates the transparency process.

The sample might be expanded 10-30% in one direction after incubation.

- 6) Perform deep imaging of the transparent brain using an appropriate dipping objective lens. Use SCALEVIEW-A2 as the immersion medium.
If necessary, immobilize the sample to the bottom of the imaging vessel with agarose.
- 7) The transparent brain can be stored for a long time in SCALEVIEW-A2 at 4 °C.

【Caution】

- 1) Ingredients may precipitate on storage at temperatures below 0 °C. Staying at room temperature (about 25 °C) returns the agent dissolved without deterioration of performance.
- 2) After the experiments, remove the residual solution from the objective lens by wiping with gauze moistened with absolute alcohol.
- 3) Be aware that the solution may cause metal corrosion (solid iron, copper, gunmetal, aluminum, etc.).
- 4) At wavelengths over 1,300 nm images may appear.

【Storage】

Store below 25 °C in the dark.

【Package】

500 mL

【Related Products】

Code No.	Description	Size
160-16061	Paraformaldehyde	100 g
164-25511	1 × PBS(-)	5 L
194-00011	Sucrose	100 g

【Reference】

- 1) Hama, H. *et al.* : *Nat Neurosci*, **14**, 1481 (2011).

FUJIFILM Wako Pure Chemical Corporation

1-2, Doshomachi 3-Chome, Chuo-Ku, Osaka 540-8605, Japan
Telephone : + 81-6-6203-3741
Facsimile : + 81-6-6201-5964
<http://www.wako-chem.co.jp>

FUJIFILM Wako Chemicals U.S.A. Corporation

1600 Bellwood Road
Richmond, VA 23237
U.S.A.
Telephone : + 1-804-271-7677
Facsimile : + 1-804-271-7791
<http://www.wakousa.com>

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