

FUJIFILM**Wako**Code No. 038-25661 (25 sheets)
034-25663 (1 roll)
032-25664 (2 sheets)**ClearTrans® PVDF Membrane,
Hydrophobic, 0.45 μm**

ClearTrans® PVDF Membrane is a versatile polyvinylidene fluoride membrane designed for blotting applications. It is more highly resistance to organic solvents and more suitable for stripping for reprobing than nitrocellulose membrane.

Product	Code No.	Size	Quantity
ClearTrans® PVDF Membrane, Hydrophobic, 0.45 μm	038-25661	10cm×10cm	25 sheets
	034-25663	30cm×400cm	1 roll
	032-25664	10cm×10cm	2 sheets

【Storage】

At room temperature

【Precautions for Use】

- Wear gloves when handling the membrane.

〈Protocol for Western Blot Transfer〉

1. Cut membrane to required dimensions.
[Note] The membrane is placed between blue protection papers. Dispose of the protection papers.
2. Wet the membrane in ethanol for 15 seconds.
3. Wash the membrane in pure water for 15 seconds.
4. Wet the membrane in transfer buffer for 5 minutes.
[Note] If discontinuous transfer buffer is used, wet the membrane in the same transfer buffer as equilibrating gel for 5 minutes.
5. Transfer proteins onto the membrane according to protocol for a blotting instrument.

FUJIFILM Wako Pure Chemical Corporation1-2, Doshomachi 3-Chome, Chuo-Ku, Osaka 540-8605, Japan
Telephone : + 81-6-6203-3741
Facsimile : + 81-6-6201-5964
<http://www.wako-chem.co.jp>**FUJIFILM Wako Chemicals U.S.A. Corporation**1600 Bellwood Road
Richmond, VA 23237
U.S.A.
Telephone : + 1-804-271-7677
Facsimile : + 1-804-271-7791
<http://www.wakousa.com>**FUJIFILM Wako Chemicals Europe GmbH**Fuggerstrasse 12
D-41468 Neuss
Germany
Telephone : + 49-2131-3111-0
Facsimile : + 49-2131-311100
<http://www.wako-chemicals.de>コード No. 038-25661 (25 枚)
034-25663 (1 巻)
032-25664 (2 枚)**クリアトランス® PVDF メンブレン,
疎水性, 0.45 μm**

クリアトランス® PVDF メンブレンは、プロットングで汎用されるポリビニリデンフルオライド性のメンブレンです。ニトロセルロースメンブレンよりも高い有機溶媒耐性を有しているため、ストリッピングバッファを用いたリプロービングにも使用できます。

品名	コード No.	大きさ	数量
クリアトランス® PVDF メンブレン, 疎水性, 0.45 μm	038-25661	10cm×10cm	25 枚
	034-25663	30cm×400cm	1 巻
	032-25664	10cm×10cm	2 枚

【保存】

室温

【ご使用前に】

- メンブレンに触れる際は手袋を着用してください。

〈使用方法〉

1. メンブレンを必要なサイズにカットしてください。
[備考] メンブレンは青色の保護紙に挟まれています。保護紙は廃棄してください。
2. カットしたメンブレンをエタノールに 15 秒浸してください。
3. カットしたメンブレンを精製水で 15 秒洗浄してください。
4. メンブレンを転写バッファに 5 分間浸してください。
[備考] 不連続バッファをご使用の場合は、ゲルの平衡化に使用する転写バッファに 5 分間浸してください。
5. プロットング装置の説明書に従って転写を行ってください。

製造発売元

富士フイルム 和光純薬株式会社
大阪市中央区道修町三丁目1番2号
Tel : 06-6203-3741

1908KA3