FUJIFILM

Wako

Code No. 016-24901 (10 g)
012-24903 (50 g)

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for Vitamin B₁ and B₂ Determination
ビタミンB₁・B₂定量用

Source : *Aspergillus oryzae*

Appearance : Slightly yellow ~ brown, powder

Potency : min. 100units/g

[Unit Definition]

One unit is that amount of enzyme needed to liberate 1 mg of phosphorus from Beta-Glycerophosphoric Acid Disodium Salt solution in 30 min. at pH 3.5 at 40 degrees C.

[General Procedure]

A. Preparation of enzyme solution

1. Add about 1g of Acid phosphatase per 100 mL pure water or buffer. Prepare just before usage.
2. Adjust pH around 4.5 and leave it at rest in cool place for a while.
3. Filter or centrifuge it and keep the supernatant for next step.

B. Digestion

Extract a sample from tissue and prepare sample including about 10 µg of Vitamin B₁ or B₂.

*For a sample from dried yeast or plant tissue like soybeans

1. Add 0.5-1 mL of 1% Acid phosphatase solution.
2. Incubate it at 40 degrees C for about 5 hours.

*For a sample from animal tissue like Porcine meat

1. Add about 10 mL of 1% Acid phosphatase solution.
2. Incubate it at 40 degrees C for about 10 hours.

Please determine specific condition for your sample, if you need accuracy in your research, because the best protocol depends on the type of sample and content of the bound form of Vitamin B₁ or B₂.

[Storage] Store at 2-10 degrees C in the dark.

[Package]

Code No.	Packaging
016-24901	10 g
012-24903	50 g

[Note] Please put the cap back or seal the bottle immediately after use, since it easily absorbs moisture.

FUJIFILM Wako Pure Chemical Corporation

1-2, Doshomachi 3-Chome, Chuo-Ku, Osaka 540-8605, Japan
Telephone : + 81-6-6203-3741
Facsimile : + 81-6-6201-5964
<http://fwk.fujifilm.co.jp>

FUJIFILM Wako Chemicals U.S.A. Corporation

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FUJIFILM Wako Chemicals Europe GmbH