FUJIFILM

Wako

Code No. 022-16231 (100 mL)

BCIP/NBT Solution

BCIP/NBT 溶液

This product contains 5-Bromo-4-chloro-3-indolylphosphate (BCIP) and Nitroblue tetrazolium (NBT).

Unreacted substrate should be colorless to light yellow in appearance. When this product is reacted with phosphatase, an insoluble permanent dark blue-purple reaction product is obtained.

This product is not recommended for microwell or immunohistochemical applications.

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[Usage]

This product is supplied as a ready-to-use solution.

This product should be allowed to equilibrate to room temperature (25℃) prior to use. In membrane assays, use enough substrate solution to completely cover the membrane surface. When the substrate system reacts with phosphatase, a blue-purple reaction product is obtained. To stop the reaction, rinse membrane in distilled water.

The reaction should be monitored and read before background color becomes too intense resulting in diminished contrast between positive and background staining.

Dilution of the substrate is not recommended. To reduce the intensity of a reaction, it is recommended that the antibody or conjugates be diluted.

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[Expiration date] Indicated on the label.

[Storage] Store at 2～10℃ in the dark.

[Package] 100 mL

* : **[Trans]** is the Japanese translation.

FUJIFILM Wako Pure Chemical Corporation

1-2, Doshomachi 3-Chome, Chuo-Ku, Osaka 540-8605, Japan
Telephone : + 81-6-6203-3741
Facsimile : + 81-6-6201-5964
<http://fwk.fujifilm.co.jp>

FUJIFILM Wako Chemicals U.S.A. Corporation

1600 Bellwood Road
Richmond, VA 23237
U.S.A.
Telephone : + 1-804-271-7677
Facsimile : + 1-804-271-7791
<http://www.wakousa.com>

FUJIFILM Wako Chemicals Europe GmbH

Fuggerstrasse 12
D-41468 Neuss
Germany
Telephone : + 49-2131-3111-0
Facsimile : + 49-2131-311100
<http://www.wako-chemicals.de>