

FUJIFILM

Wako

Code No. 045-35121 (1 mg)
041-35123 (5 mg)
049-35124 (25 mg)

3 × DYKDDDDK Peptide 3 × DYKDDDDK ペプチド

The 3×DYKDDDDK Peptide is used for the competitive elution of MDYKDHG DYKDHDI DYKDDDDK (3×DYKDDDDK) tag or DYKDDDDK tag fusion proteins from the immune complex after immunoprecipitation by using anti 3×DYKDDDDK or DYKDDDDK tag antibody.

This product is for laboratory use only : use in any such application is the responsibility of the user.

【Formulation】

Trifluoroacetic acid salt.

【Amino acid sequence】

N-Met-Asp-Tyr-Lys-Asp-His-Asp-Gly-Asp-Tyr-Lys-Asp-His-Asp-Ile-Asp-Tyr-Lys-Asp-Asp-Asp-Lys-C

【Preparation of stock solution】

Dissolve in 1×TBS (50 mmol/L Tris-HCl, pH 7.4, 150 mmol/L Sodium Chloride) to a final concentration of 1 mg/mL. Aliquot and store at −20°C. Avoid repeated freeze and thaw cycle.

【Working concentration】

Standard : 100 μg/mL in 1×TBS

A working concentration depends on the experiment.

【Storage】

Store at −20°C and protect from light.

【Package】

Code No.	Packaging
045-35121	1 mg
041-35123	5 mg
049-35124	25 mg

FUJIFILM Wako Pure Chemical Corporation

1-2, Doshomachi 3-Chome, Chuo-Ku, Osaka 540-8605, Japan
Telephone : + 81-6-6203-3741
Facsimile : + 81-6-6201-5964
<http://fwk.fujifilm.co.jp>

Group Companies



Distributors



2407KA1