

**FUJIFILM**

**Wako**

Code No. 043-31381 (1 g)  
049-31383 (5 g)

## Deoxyribonucleic Acid, from Salmon Sperm

### デオキシリボ核酸, サケ精液由来

This product is the sheared DNA from salmon sperm. It is used to prevent non-specific binding of the probe to the membrane in Southern and Northern Hybridization. This product should be sonicated, extracted with phenol/chloroform and precipitated with ethanol before using.

#### 【Appearance】

Lyophilized

#### 【Size of DNA】

Approximately 300 bp ~ 1,000 bp

#### 【Storage】

2 ~ 10°C・遮光保存

#### 【Package】

Code No.	Package
043-31381	1 g
049-31383	5 g

## FUJIFILM Wako Pure Chemical Corporation

1-2, Doshomachi 3-Chome, Chuo-Ku, Osaka 540-8605, Japan  
Telephone : + 81-6-6203-3741  
Facsimile : + 81-6-6201-5964  
<http://fwk.fujifilm.co.jp>

Group Companies



Distributors

