

FUJIFILM

Wako

Code No. 088-10531 (1 mg)
084-10533 (5 mg)

6×Histidine Peptide 6×ヒスチジンペプチド

The 6×Histidine (His) Peptide is synthesized artificially and consisted of 6 amino acids. The 6×His Peptide is used for the competitive elution of 6×His tag fusion proteins from anti 6×His antibody.

This product is for laboratory use only ; use in any such application is the responsibility of the user.

【Formulation】

Trifluoroacetic acid salt.

【Amino acid sequence】

His-His-His-His-His-His

【Preparation of stock solution】

Dissolve in 1×TBS (10 mmol/L Tris-HCl, pH 7.4, 150 mmol/L Sodium Chloride) to a final concentration of 5-10 mg/mL. Aliquot and store at -20°C. Avoid repeated freeze and thaw cycle.

【Working concentration】

Standard : 100-150 µg/mL in 1×TBS

A working concentration depends on the experiment.

【Storage】

Store at -20°C and protect from light.

Keep container tightly closed in a dry and well-ventilated place.

【Package】

Code No.	Packaging
088-10531	1 mg
084-10533	5 mg

FUJIFILM Wako Pure Chemical Corporation

1-2, Doshomachi 3-Chome, Chuo-Ku, Osaka 540-8605, Japan
Telephone : + 81-6-6203-3741
Facsimile : + 81-6-6201-5964
<http://fwk.fujifilm.co.jp>

FUJIFILM Wako Chemicals U.S.A. Corporation

1600 Bellwood Road
Richmond, VA 23237
U.S.A.
Telephone : + 1-804-271-7677
Facsimile : + 1-804-271-7791
<http://www.wakousa.com>

FUJIFILM Wako Chemicals Europe GmbH

Fuggerstrasse 12
D-41468 Neuss
Germany
Telephone : + 49-2131-311-0
Facsimile : + 49-2131-311100
<http://www.wako-chemicals.de>

2306KA2