FUJIFILM**Wako**

For Research Use Only. Not for use in diagnostic procedures.

Code No. 122-07031 (80 μ L)**LBIS™ Porcine Insulin Standard Solution****Volume**80 μ L/bottle**Insulin Concentration**

240 ng/mL

Purification Method

Reversed-phase chromatography

Buffer

Phosphate buffer (with stabilizer)

Storage

2°C - 10°C

Procedure

Porcine insulin can be quantitatively measured by replacing the Insulin Standard included in the LBIS™ Rat Insulin ELISA Kit (Code No. 299-94801) with this product. When using a portion of this product, remove it from the refrigerator immediately before use to prepare the dilution. Immediately and tightly close the remaining stock solution without allowing it to return to room temperature and store it at 2°C - 10°C. Use the prepared dilutions promptly and do not store them.

Porcine Insulin Standard Solution (240 ng/mL) for Standard Curve Preparation

Prepare standard solutions using the LBIS™ Porcine Insulin Standard Solution (240 ng/mL, stock solution) and (C) Buffer supplied with the kit. The following is an example. *Conversion to μ IU/mL is calculated using 26 IU/mg.

Standard Solution Volume	(C) Buffer	Concentration (ng/mL)	Concentration (μ IU/mL*)
10 μ L of Standard stock solution	190 μ L	12	312
100 μ L of 12 ng/mL solution	100 μ L	6.0	156
100 μ L of 6.0 ng/mL solution	100 μ L	3.0	78
100 μ L of 3.0 ng/mL solution	100 μ L	1.5	39
100 μ L of 1.5 ng/mL solution	100 μ L	0.75	19.5
100 μ L of 0.75 ng/mL solution	100 μ L	0.375	9.75
100 μ L of 0.375 ng/mL solution	100 μ L	0.188	4.88
0 (Blank)	100 μ L	0	0

Expiration Date

6 months from the date of manufacture (when unopened and kept refrigerated). The expiration date is indicated on the outer label of the kit.

Precaution

As the Standard Solution may be dispersed inside the bottle, collect it at the bottom by centrifugation or similar methods before opening the cap.

(Recommended centrifugation conditions : 500 g - 1,000 g, 10 seconds - 15 seconds, 2°C - 10°C)

LBIS™ Porcine Insulin Standard Solution

[Storage] Store at 2°C - 10°C

[Expiration Date] Indicated on the label.

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