FUJIFILM**Wako**

<for Research Use Only> Code No. 126-06711 (100 tests)

Leucomalachite Green Reaction Reagent [S Leuco]

Leucomalachite Green is the reduced form of Malachite Green, and is converted from colorless to blue-green (Malachite Green) by oxidation.

This oxidation reaction is catalyzed by hemoglobin under an acidic condition to produce blue-green color, and often used to detect the presence of faint blood stains.

General Leucomalachite Green reaction reagents need to be reconstituted at the time of use, because prepared solution is less stable, and cannot be stored for a long time after preparation.

This Ready-to-Use product contains all the necessary reagents premixed for Leucomalachite Green reaction, and does not need solution preparation.

【Additional required reagent and material】

Reagent :

- 1 mol/L Citric acid solution
ex) Prepare 100 mL of 1 mol/L Citric acid solution.
Dissolve 19.21 g of Citric acid or 21.01 g of Citric acid monohydrate with distilled water. Make up to 100 mL.

Material :

- Cotton swab or equivalents

【Usage】

1. Apply citric acid solution to the tip of the cotton swab.
2. Wipe a specimen with a cotton swab moistened with citric acid solution.
3. Push the cotton swab with which specimen was wiped away to this reagent.
4. Check the color development of this reagent.

【Precaution for Use】

- Wear safety glasses, protective gloves, and protective mask when used.
- Please apply citric acid solution to the tip of cotton swab immediately before specimen collection.
- Room temperature is defined as 30 °C or less. Avoid storing under an environment that is excessively hot.
- Wipe specimen with the tip of the cotton swab to enable steady contact between the specimen and this reagent.

- Since color development of this reaction disappears after a long period of time, check the presence or absence of coloring immediately after contacting the specimen with this reagent.

【Storage】

Store at room temperature. Protect from light.

【Package】

100 tests

【Reference】

Cox, M., *J. Forensic Sci.*, **36** (5), 1503 (1991).

FUJIFILM Wako Pure Chemical Corporation

1-2, Doshomachi 3-Chome, Chuo-Ku, Osaka 540-8605, Japan
Telephone : + 81-6-6203-3741
Facsimile : + 81-6-6201-5964
<http://fwk.fujifilm.co.jp>

FUJIFILM Wako Chemicals U.S.A. Corporation

1600 Bellwood Road
Richmond, VA 23237
U.S.A.
Telephone : + 1-804-271-7677
Facsimile : + 1-804-271-7791
<http://www.wakousa.com>

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