LBIS™ Mouse Albumin Control Serum

Please make sure you read this manual before using the product.

【Intended Use】

Use this product as a positive control for the LBIS™ Mouse Urinary Albumin Assay Kit.

【Contents】

LBIS™ Mouse Albumin Control Serum, 1 mL/bottle

(The concentration varies by lot.)

【Procedure】

Thaw immediately before the assay and use as a positive control for the LBIS™ Mouse Urinary Albumin Assay Kit (Code No. 295-92701/291-92703) without dilution. After thawing, store at 2°C - 10°C and use within 7 days. Avoid repeated freeze-thaw cycles.

【Precautions/Restrictions】

1. This product is a positive control specifically designed for FUJIFILM Wako kits. Do not use it with kits or reagents from other manufacturers.
2. Store this product under the specified conditions. Do not use products that have passed their expiration date.
3. Do not use this product for any purposes other than those described in this manual.
4. The concentration listed on the label varies by lot.
5. Avoid repeated freeze-thaw cycles.

【Safety Precautions】

1. If the product comes into contact with the human body, immediately rinse the affected area with plenty of water and seek medical attention as needed.
2. Since it contains serum components, when handling this product, take the same precautions that apply to clinical specimens due to the potential risk of viral infection.
3. Dispose of the product in accordance with applicable laws and regulations related to waste management and wastewater standards.

LBIS™ Mouse Albumin Control Serum

[Storage]	Store frozen (-20°C or below), protect from light.
[Indicated Value]	Indicated on the label.
[Expiration Date]	Indicated on the label.
[Lot #]	Indicated on the label.
[Package]	1 mL × 10 bottles.
[Cat #]	127-07081

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