LabAssay™ GLDH Control

Please, read this instruction carefully before use.

【Intended Use】

This product is used as a control for LabAssay™ GLDH (AA-type).
(Code No. 291-97801).

【Volume】

LabAssay™ GLDH Control 500 μ L/bottle
(The displayed value of GLDH differs for each lot.)

【How to Use】

Allow the reagent to thaw completely at room temperature before use. Use the solution undiluted as the detection sample exclusively for LabAssay™ GLDH (AA-type).

After thawing, gently mix by inversion to avoid foam formation; use of a rolling mixer or gentle pipetting is recommended.

Avoid vigorous agitation, as it may result in decreased enzyme activity.

Use the reagent on the day of thawing. Repeated freeze-thaw cycles are strictly prohibited.

When measuring LabAssay™ GLDH Control as a sample, the expected measurement value falls within $\pm 15\%$ of the stated value.

【Important Notice in the Treatments】

1. This product is a control for use with LabAssay™ GLDH (AA-type) only. Do not use it with other companies' kits or reagents.
2. Please store this product under the specified conditions and do not use it after the expiration date.
3. Do not use this product for any purpose other than those described in this manual.
4. The displayed value of GLDH differs for each lot.
5. Do not repeat freezing and thawing.

【Precaution】

1. Avoid contact with skin. In case of contact, wash it thoroughly with water and seek medical attention if needed.
2. This product is derived from serum, therefore, handle this product as potentially bio-hazardous and treat it carefully as your samples.
3. Dispose used and unused controls in accordance with applicable national or regional regulatory rules for discard of biological wastes.

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