

FUJIFILM

Wako

Code No. 193-16375 (500 mL)

Stripping Solution

Stripping Solution is used to remove primary and secondary antibodies from a Western blot membrane. Stripping is useful for investigating more than one protein on the same membrane, to save time and conserve samples.

Protocol

1. Wash the already probed membrane with TBS-T or PBS-T for 5 minutes two times.
2. Incubate the membrane in Stripping Solution with agitation for 10 minutes at room temperature. Use a sufficient volume that will cover the membrane.
3. Remove the membrane from the Stripping Solution and wash three times with TBS-T or PBS-T for 5 minutes.
4. Perform antigen-antibody reaction again.

Notes

- Stripping Solution should not be used with tweezers and a tray made of metal.
- Use Stripping Solution with membranes detected by chemiluminescent detection reagents. It is not recommended to use Stripping Solution with chromogenic detection reagents because it leaves a visible stain on the membrane and Stripping Solution cannot strip it.

Package : 500 mL

Storage : at room temperature

FUJIFILM Wako Pure Chemical Corporation

1-2, Doshomachi 3-Chome, Chuo-Ku, Osaka 540-8605, Japan
Telephone : + 81-6-6203-3741
Facsimile : + 81-6-6201-5964
<http://fwk.fujifilm.co.jp>

FUJIFILM Wako Chemicals U.S.A. Corporation

1600 Bellwood Road
Richmond, VA 23237
U.S.A.
Telephone : + 1-804-271-7677
Facsimile : + 1-804-271-7791
<http://www.wakousa.com>

FUJIFILM Wako Chemicals Europe GmbH

Fuggerstrasse 12
D-41468 Neuss
Germany
Telephone : + 49-2131-3111-0
Facsimile : + 49-2131-311100
<http://www.wako-chemicals.de>

コード No. 193-16375 (500 mL)

ストリッピング溶液

ストリッピング溶液は、ウエスタンブロットを行ったメンブレンから抗体を除去する試薬です。同一メンブレン上で複数の抗原検出を可能にし、時間短縮およびサンプルを節約できます。

使用方法

1. 抗原抗体反応を行ったメンブレンをTBS-TかPBS-Tで5分間洗浄します。2回洗浄します。
2. ストリッピング溶液にメンブレンを浸しながら、室温で10分間振とうします。十分にメンブレンが浸る量のストリッピング溶液を使用します。
3. ストリッピング溶液からメンブレンを取り出し、TBS-TかPBS-Tで5分間洗浄します。3回洗浄します。
4. 再度、抗原抗体反応を行ってください。

備 考

- 金属製のピンセットやトレイはなるべく使用しないでください。
- 化学発光試薬で検出したメンブレンに使用してください。発色基質で検出すると、色素がメンブレンに沈着し、本品で剥がすことはできません。

容 量 : 500 mL

保存条件 : 室温

製造発売元

富士フイルム 和光純薬株式会社
大阪市中央区道修町三丁目1番2号
Tel : 06-6203-3741