FUJIFILM

Wako

Code No. 193-19533 (10 mL × 100)
191-19534 (25 mL × 50)
197-19536 (30 mL × 50)


SafeCap™ Wako 10% Formalin Neutral Buffer Solution

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Tissue fixation is an important procedure to protect biological specimens from degradation due to autolysis and decay, ensuring stable preservation of proteins and other components within the tissues. Formalin fixation is widely used in research and pathology, including microscopy specimen preparation. SafeCap™ Wako is a container that allows tissue samples to be safely fixed and preserved without exposure to formalin.

【Features】

- The cap contains formalin solution, so it is ready to use immediately.
- Easy operation: simply place the sample in the container and tighten the cap. Formalin solution can be safely filled into the container.
- A wide-mouth container that makes it easy to put in samples.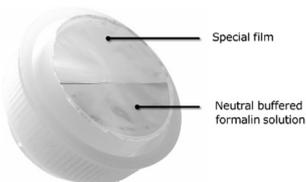

【Storage】

Store at room temperature in the dark.

【Procedure】

① Open the cap.

② Place the sample in the container.

③ Tighten the cap until it cannot be turned any further.

④ Slowly tilt the container and check that the neutral buffered formalin solution flows into the container.

*The liquid flows due to air exchange inside the container and cap. Do not shake vigorously or tighten the cap excessively, as this may cause breakage or leakage.

【Precautions】

- Avoid direct sunlight and store in a dark place.
- When handling, wear a gas mask or hose mask, protective gloves, and glasses as necessary.
(For details, please refer to the handling precautions on the label.)

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