FUJIFILM**Wako**

Code No. 199-18153 (1 mL)

For Genetic Research

SAFELOOK™ Load-Green (6×)

SAFELOOK™ Load-Green (6×) is an non-carcinogenic and safe nucleic acid stain for the visualization of double-strand DNA, single-strand DNA, and RNA in agarose and polyacrylamide gels. This product is developed to replace toxic Ethidium Bromide, and the Ames-test was conducted to confirm its safety. SAFELOOK™ Load-Green (6×) is provided in a form of 6× sample loading dye and is to be added to your samples only.^{※1}

※1 Please optimize volume of reagent, amount of nucleic acid, thickness of the gel, and staining time if necessary.

【Features】

- ・ Substitute for 6× sample loading dye
- ・ Dye emits green fluorescence
- ・ Visualize dsDNA, ssDNA, and RNA
- ・ Excitation spectra (490 nm)
- ・ Emission spectra (525 nm)
- ・ Sensitivity range is between 0.2-0.6 ng
- ・ Work with UV and LED transilluminator
- ・ Available for real-time visualization

【Protocol】

- 1) Prepare agarose gels or polyacrylamide gels.
- 2) Mix samples and DNA marker with SAFELOOK™ Load-Green (6×) at a 1 : 5 (dye : sample) dilution rate.
- 3) Following electrophoresis, view the results under UV or LED light.

【Storage】

2-10℃

【Package】

Code No.	Packaging
199-18153	1 mL

FUJIFILM Wako Pure Chemical Corporation

1-2, Doshomachi 3-Chome, Chuo-Ku, Osaka 540-8605, Japan
Telephone : + 81-6-6203-3741
Facsimile : + 81-6-6201-5964
<http://fwk.fujifilm.co.jp>

FUJIFILM Wako Chemicals U.S.A. Corporation

1600 Bellwood Road
Richmond, VA 23237
U.S.A.
Telephone : +1-804-271-7677
Facsimile : +1-804-271-7791
<http://www.wakousa.com>

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