

FUJIFILM

Wako

Code No. 201-21271 (1 mL)

TamavidinTM2-REV, recombinant, Solution タマビジンTM2-REV, 組換え体, 溶液

TamavidinTM2 is an avidin-like protein cloned from *Pleurotus cornucopiae*, a species of mushroom. By introducing a point mutation into the domain of TamavidinTM2 involved in binding to biotin, a new protein was created that can reversibly bind to biotin. This protein was named TamavidinTM2-REV.

【Features】

- Able to reversibly bind to biotin
: The biotinylated factor can be recovered from TamavidinTM2-REV by adding excess biotin.
- Near-neutral isoelectric point (pI)
: Non-specific binding due to charge is prevented.
- Protease-resistant
: It can be treated with protease after it has bound the biotinylated factor.

【Source】

E. coli expressed mushroom (*Pleurotus cornucopiae*) TamavidinTM2-REV.

【Product concentration】

Approximately 1 mg/mL

【Formulation】

20 mmol/L potassium phosphate buffer (pH 7.0)

【Storage conditions】

Store at -20°C.

【Molecular weight】

Approximately 60 k (SDS-PAGE) (tetramer)

【Isoelectric point (pI)】

Approximately 7.2

【Thermal stability】

*T*_m = Approximately 80°C

【An example of usage】 [Recovery of biotinylated factor]

Immobilize TamavidinTM2-REV on the bead carrier (e.g., agarose, magnetic beads)



Mix it with the biotinylated factor and react



Remove the supernatant by centrifugation
or using a magnetic stand
↓ Add washing buffer
↓ Mix
Remove the supernatant by centrifugation
or using a magnetic stand
↓ Add biotin solution* as an elution buffer
↓ Mix
Recover the supernatant as a purified fraction by centrifugation
or using a magnetic stand

} Repeat multiple times

* Examples of how to prepare biotin solution as an elution buffer

- 2 mmol/L biotin solution
Use PBS to make a biotin concentration of 2 mmol/L.
- 20 mmol/L biotin solution
Use 50 mmol/L potassium phosphate buffer (pH 7.0) to make a biotin concentration of 20 mmol/L.

【References】

- 1) Takakura, Y. *et al.*, "Tamavidin 2-REV : An engineered tamavidin with reversible biotin-binding capability.", *J. Biotechnol.*, **164**(1), 19-25 (2013).
- 2) Motani, K. *et al.*, "BioID screening of biotinylation sites using the avidin-like protein Tamavidin 2-REV identifies global interactors of stimulator of interferon genes (STING).", *J. Biol. Chem.*, **295**(32), 11174-11183 (2020).

【Package】

1 mL

FUJIFILM Wako Pure Chemical Corporation

1-2, Doshomachi 3-Chome, Chuo-Ku, Osaka 540-8605, Japan
Telephone : + 81-6-6203-3741
Facsimile : + 81-6-6201-5964
<http://ffwk.fujifilm.co.jp>

FUJIFILM Wako Chemicals U.S.A. Corporation

1600 Bellwood Road
Richmond, VA 23237
U.S.A.
Telephone : + 1-804-271-7677
Facsimile : + 1-804-271-7791
<http://www.wakousa.com>

FUJIFILM Wako Chemicals Europe GmbH

Fuggerstrasse 12
D-41468 Neuss
Germany
Telephone : + 49-2131-3111-0
Facsimile : + 49-2131-311100
<http://www.wako-chemicals.de>