LabAssay™ Albumin

Please read this instruction carefully before use.

1. Intended use

This product is a research-use reagent for the measurement of albumin using the bromocresol green (BCG) method.

Albumin is a plasma protein synthesized in the liver. Serum or plasma albumin levels are important indicators in the diagnosis and management of liver disease and hepatic dysfunction. In addition, kidney diseases characterized by protein leakage into the urine are known to cause decreases in blood total protein and albumin levels, leading to hypoalbuminemia.

The albumin/globulin ratio (A/G ratio) can be calculated from the measured albumin and total protein concentrations and serves as an adjunct diagnostic indicator for liver disease, kidney disease, nutritional disorders, multiple myeloma, and chronic inflammatory diseases. Albumin and total protein are commonly measured together.

This product includes albumin standards derived from bovine, mouse, and rat. Albumin measurement using the BCG method is known to exhibit species-dependent differences in reactivity. Using the bovine, mouse, and rat albumin standards provided in this kit enables more accurate quantification of albumin concentrations.

2. Storage and expiration date

Store at 2°C - 10°C and do not freeze.

The expiration date is indicated on the label on the outer box of the kit.

3. Kit Component Reagents

	Components	Use Status	Amount
(1)	Chromogen Reagent	Solution	65 mL/1 bottle
(2)	Bovine Albumin Standard	Freeze-dried	1 bottle
(3)	Mouse Albumin Standard	Freeze-dried	1 bottle
(4)	Rat Albumin Standard	Freeze-dried	1 bottle

* For 250 tests

4. Principle of the method

When the Chromogen Reagent is added to a sample, albumin binds to bromocresol green (BCG) to form a blue-colored complex. Albumin concentration is determined by measuring the intensity of the blue color.

5. Equipment or supplies required but not provided in the kit

- 96-well microplate (transparent type)
- Micropipette
- Microtube
- Pipette
- Plate mixer
- Microplate reader with 630 nm wavelength filter

6. Preparation of reagents

① Chromogen Reagent : Ready to use.

② Bovine Albumin Standard : Add the specified volume* of purified water as indicated in the "Reconstitution of standard" section to the standard vial. Allow it to stand for 10 minutes, then gently mix until completely dissolved, avoiding foaming, to prepare the standard stock solution (5.00 g/dL). Thereafter, dilute with purified water at room temperature.

Store any remaining standard stock solution in aliquots frozen. It is recommended to store frozen at -35°C or below. Do not repeat freeze-and-thaw cycles.

*Refer to the "Reconstitution of standard" section on our product webpage. The volume of purified water to be added may vary depending on the lot; therefore, confirm the appropriate volume for each lot.

*Use purified water equilibrated to room temperature to dissolve the standard.

*Standard solutions diluted to each concentration should be used immediately and must not be stored.

③ Mouse Albumin Standard : For mouse samples, use the mouse albumin standard. See Section ② for preparation, dilution, and storage instructions.

④ Rat Albumin Standard : For rat samples, use the rat albumin standard. See Section ② for preparation, dilution, and storage instructions.

Example of preparation of dilution series of standard solution

Standard Concentration (g/dL)	7.5	5.0	2.5	1.25	0.625	0.0
Standard Solution (μL)	Original standard solution : 50	Original standard solution : 50	Original standard solution : 50	50 *	50 *	—
Water (μL)	—	—	50	50	50	50
Well Volume (μL)	3**	2	2	2	2	2

* : One rank higher standard solution

** : The standard solution is typically dispensed at 2 μL per well ; however, 3 μL is used for the 7.5 g/dL standard. Accordingly, the displayed value has been corrected to account for the increased volume.

7. Preparation of specimen

Serum/Plasma

- Analyze specimens immediately after collection.
- The anticoagulant heparin, EDTA · 2Na and citrate does not significantly influence the assay when used in normal amounts.
- Serum samples with a high degree of hemolysis may affect the measurement results. In such cases, collecting the blood sample again is recommended or the results should be used for reference only.
- Chyle and ascorbic acid affect the measured values.
- High levels of bilirubin in the specimen may affect the measured values.
- If the measured value exceeds the measurable range, dilute the sample with saline solution before measurement. Multiply the result by the dilution factor.

8. Assay procedure

Ensure that the reagents are brought up to room temperature (20°C - 25°C) before use.

- (1) Dispense 250 μL of Substrate-Enzyme Solution into each well of a 96-well microplate.
 - (2) Dispense 2 μL of each concentration of standard solution into the well in which the standard is to be measured. (Dispense the 7.5 g/dL standard at the volume specified in the standard dilution example.)
 - (3) Dispense 2 μL of sample into the well in which the sample is to be measured.
 - (4) Shake the 96-well microplate on a plate mixer (* ①).
 - (5) Allow to react at room temperature at 20 - 25°C for 20 minutes.
 - (6) Immediately after reaction, measure the absorbance of each well at 630 nm using a spectrophotometer for microplates.
- (* ①) Guideline for shaking : 600 rpm - 700 rpm for 10 seconds, repeated 3 times.

9. Calculation


- (1) Create a calibration curve for each measurement, with the standard concentration (g/dL) on the X-axis and absorbance on the Y-axis.
- (2) From the calibration curve, read the albumin (g/dL) corresponding to the absorbance of the sample. If a diluted sample is used, the concentration reading is multiplied by the sample dilution ratio to obtain the measurement value.

* If the absorbance of a sample is outside the standard curve, dilute the sample with saline to an appropriate dilution ratio and repeat the assay.

* In computer software calculations, we recommend using 4 parameters or 5 parameters.

10. Standard Curve (example)

- Calibration range : 0.625 ~ 7.5 g/dL

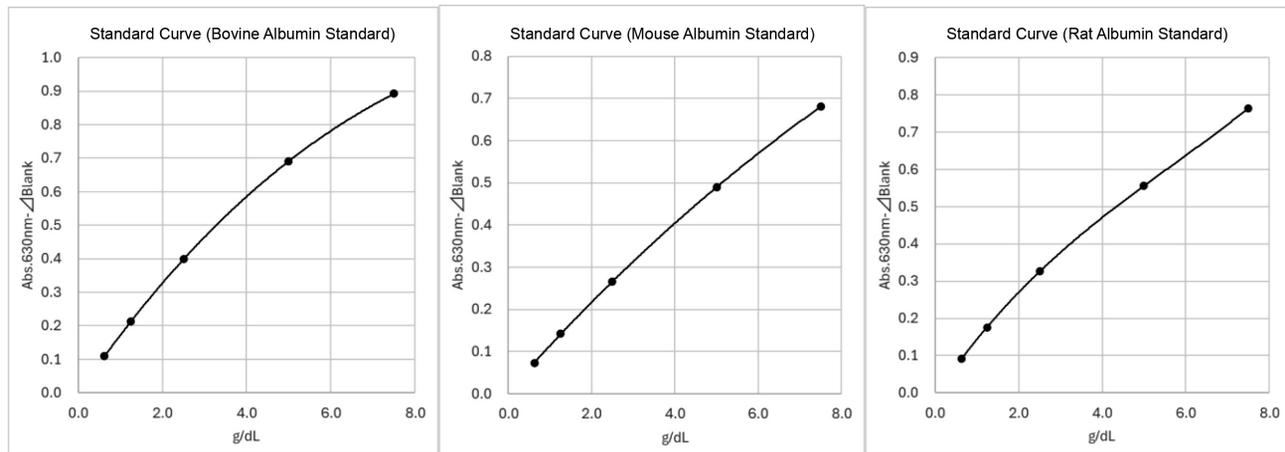11. Performance

- Accuracy
- When control serum of known concentration is measured, the measured value is within $\pm 15.0\%$ of the known concentration.
- Repeatability
- When the same sample is measured simultaneously 5 times, the CV (%) of measured values is 15.0% or less.

Notes

- Avoid using this product at temperatures or reaction times other than those specified.
- The color development is stable, and the absorbance shows almost no change even after standing at room temperature for 1 hour.
- If any reagent accidentally comes into contact with the mouth, eyes, or skin, flush immediately with a large amount of water. Consult a physician if necessary.
- When using a pipette, do not aspirate by mouth; always use a safety pipette filler or equivalent device.
- Store reagents under the specified conditions and do not use expired reagents.
- After opening, use the reagent as soon as possible. If it is stored, close the lid and store under the specified conditions.
- Do not use the container or accessories included with this product for anything other than their intended purposes.
- Handle samples being cognizant of the risk of infection (e.g., viruses).
- Wear disposable gloves to avoid the risk of infection when performing a test.
- Be careful regarding contamination when collecting reagents with a pipette.
- Do not mix and use reagents with different lot numbers.
- When disposing of the reagents, dispose of them according to Waste Management and Public Cleansing Law (Waste Disposal and Cleaning Act) and wastewater standards. The chromogenic solution contains bromocresol green, a phenol compound, at 120 mg/L.
- Any reagent or reagent bottle that comes into contact with a specimen should be considered potentially infectious and handled accordingly.
- Any scattering/splashes of samples or waste fluids should be removed with a disinfectant such as sodium hypochlorite (effective chlorine concentration 1,000 ppm)

LabAssay™ Albumin

[Storage] Store at 2°C - 10°C (Do not freeze)

[Term of validity] Indicated on the label.

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