## EV-Perm™ Permeabilization Pretreatment Kit for Exosome Membrane

Please visit our website through URL below and enter product code (294-85701) to learn more about product features and instructions.

<https://labchem-wako.fujifilm.com>

### 【Product Information】

| Code No.  | Product Name                                                    | Volume | Storage       |
|-----------|-----------------------------------------------------------------|--------|---------------|
| 294-85701 | EV-Perm™ Permeabilization Pretreatment Kit for Exosome Membrane | 1 kit  | Frozen (-20℃) |

### 【Kit Contents】

| Reagent          | Volume   | Storage                                                             |
|------------------|----------|---------------------------------------------------------------------|
| Reagent A (20×)  | 1×500 μL | Frozen (-20℃)                                                       |
| Reagent B (10×)  | 1×5 mL   | Frozen (-20℃) before thawing<br>Refrigeration (2–10℃) after thawing |
| Reagent C (100×) | 1×500 μL | Frozen (-20℃) before thawing<br>Refrigeration (2–10℃) after thawing |

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EV-Perm™ Permeabilization Pretreatment Kit for Exosome Membrane (EV-Perm™) enables the detection of internal extracellular vesicle proteins when PS Capture™ Exosome ELISA Kit series or PS Capture™ Exosome Flow Cytometry Kit are used in combination with EV-Perm™.

This product is research use only.

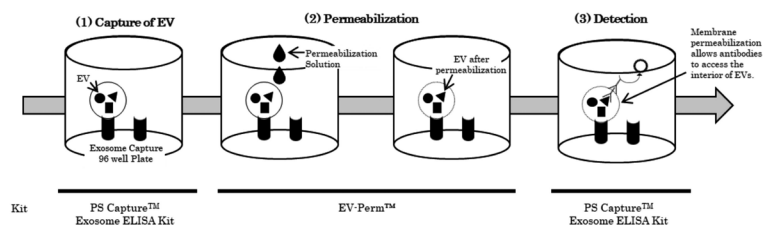
### 【Attention】

When using EV-Perm™, it is necessary to combine it with one of the following.


| Code No.  | Product Name                                       |
|-----------|----------------------------------------------------|
| 297-79201 | PS Capture™ Exosome ELISA Kit (Anti Mouse IgG POD) |
| 298-80601 | PS Capture™ Exosome ELISA Kit (Streptavidin HRP)   |
| 297-79701 | PS Capture™ Exosome Flow Cytometry Kit             |

### [Procedure for detection using EV-Perm™]

■ If you use EV-Perm™ in conjunction with the PS Capture™ Exosome ELISA Kit series

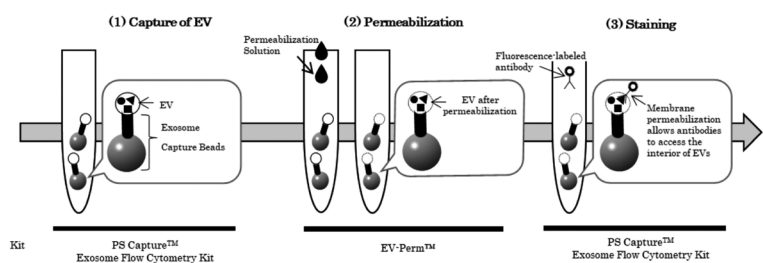■ If you use EV-Perm™ in conjunction with the PS Capture™ Flow Cytometry Kit

## FUJIFILM Wako Pure Chemical Corporation

1-2, Doshomachi 3-Chome, Chuo-Ku, Osaka 540-8605, Japan  
 Telephone : + 81-6-6203-3741  
 Facsimile : + 81-6-6201-5964  
<http://ffwk.fujifilm.co.jp>

### FUJIFILM Wako Chemicals U.S.A. Corporation

1600 Bellwood Road  
 Richmond, VA 23237  
 U.S.A.  
 Telephone : + 1-804-271-7677  
 Facsimile : + 1-804-271-7791  
<http://www.wakousa.com>

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### FUJIFILM Wako Chemicals Europe GmbH