

**LBIS™ Human Apo B-48 Control Set**

Please, read this instruction carefully before use.

LBIS™ Human Apo B-48 Control Set includes quality control sera for LBIS™ Human Apo B-48 ELISA Kit.

**【Components】**

Reagents	Use Status	Amounts		
		10 bottles × 2 sets	25 bottles × 2 sets	50 bottles × 2 sets
1) Human Apo B-48 Control-L	Frozen. Use after thawing.	250 $\mu$ L/10 bottles	250 $\mu$ L/25 bottles	250 $\mu$ L/50 bottles
2) Human Apo B-48 Control-H	Frozen. Use after thawing.	250 $\mu$ L/10 bottles	250 $\mu$ L/25 bottles	250 $\mu$ L/50 bottles

\*The concentration is indicated on the label for each lot.

**【How to use】**

Take out from freezer and thaw. Use the solutions as positive control sera as they are for LBIS™ Human Apo B-48 ELISA Kit. After thawing, keep them at 2°C - 10°C and use them up within 8 hours. Do not repeat freezing and thawing.

**【Important Notice in the Treatments】**

- (1) This LBIS™ Human Apo B-48 Control Set is intended to use for only LBIS™ Human Apo B-48 ELISA Kit. Do not use for other maker's products.
- (2) Store this product under the condition indicated below. Do not use after the expiration date.
- (3) Do not use this product for other purposes than indicated in this manual.
- (4) The concentrations of the calibrated substances will be different in each lot.
- (5) Do not repeat freezing and thawing.

**【Statements and Precaution】**

- (1) Avoid contact with skin. In case of contact, wash it thoroughly with water and seek medical attention if needed.
- (2) This product is derived from human serum, therefore, handle this product as potentially bio-hazardous and treat it carefully as your samples.
- (3) Dispose used and unused controls in accordance with applicable national or regional regulatory rules for discard of biological wastes.

**【Storage Condition】**

Store this original product in a dark freezer at below -20°C.

**【Expiration date】**

12 months from production. Expiration date is indicated on the label.

**【Unit of Package】**

LBIS™ Human Apo B-48 Control Set :

Human Apo B-48 Control-L, 250  $\mu$ L/1 bottleHuman Apo B-48 Control-H, 250  $\mu$ L/1 bottle

LBIS™ Human Apo B-48 Control Set

[Storage] Store at below -20°C.

[Expiration date] Indicated on the label

[Package] 10 bottles × 2 sets

25 bottles × 2 sets

50 bottles × 2 sets

[Cat #] 292-89009 (10 bottles × 2 sets)

298-89001 (25 bottles × 2 sets)

294-89003 (50 bottles × 2 sets)

---

## FUJIFILM Wako Pure Chemical Corporation

1-2, Doshomachi 3-Chome, Chuo-Ku, Osaka 540-8605, Japan

Telephone : + 81-6-6203-3741

Facsimile : + 81-6-6201-5964

<http://fwk.fujifilm.co.jp>

### FUJIFILM Wako Chemicals U.S.A. Corporation

1600 Bellwood Road

Richmond, VA 23237

U.S.A.

Telephone : + 1-804-271-7677

Facsimile : + 1-804-271-7791

<http://www.wakousa.com>

### FUJIFILM Wako Chemicals Europe GmbH

Fuggerstrasse 12

D-41468 Neuss

Germany

Telephone : + 49-2131-3111-0

Facsimile : + 49-2131-311100

<http://www.wako-chemicals.de>