

All Made in Japan, High-Quality JQ Vial Kit

- Japan Quality -






- Manufactured in Japan
- Low sample adsorption
- Minimized Na⁺/impurities elution

All made in Japan, High-Quality vial kit JQ Vial Kit is an "all made in Japan" vial kit comprising a glass vial and a septum cap, both made in Japan. The vial is manufactured by a unique method with an optimized processing temperature in the high-temperature molding process, enabling low adsorption and low alkaline elution without chemical treatment with silane, etc. The septum cap is also of high quality, minimizing elution of impurities.

In chromatographic analysis such as GC and LC, there are risks of unexpected analytical defects such as sample adsorption, decomposition, and appearance of ghost peaks, depending on the quality of the sample containers, the vial and septum cap. As the recent developments in analytical equipment and technology bring a need for ever more sophisticated analysis, it is critical to focus on the quality of the sample containers as well.

(1) Comparison of Sample Adsorption to Vials


Sample adsorption to vials was tested to compare our Japan-made glass vial with vial products from other companies. Diquat, used as a herbicide and known to be highly adsorptive to glass, was used as sample. After sealing diquat solution into a vial, recovery up to 12 hours was measured by HPLC.

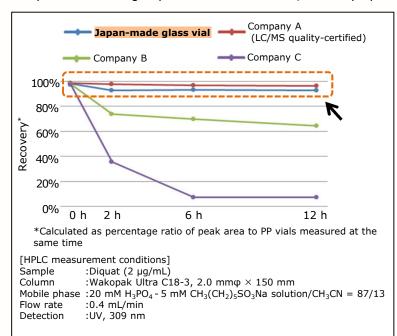

Diquat adsorption to Japanmade glass vial was hardly observed up to 12 hours, demonstrating as low adsorption as LC/MS certified high-quality glass vials from other companies. As sample adsorption to vials causes sample loss, use of lowadsorption vials is important to avoid problems in analysis.

(2) Comparison of Na⁺ Elution from Vials


Na⁺ elution from vials was tested to compare our Japan-made glass vial with vial products from other companies. After sealing purified water into vials, Na⁺, known to be an alkaline component eluted from glass, was measured by atomic absorption spectrometry.

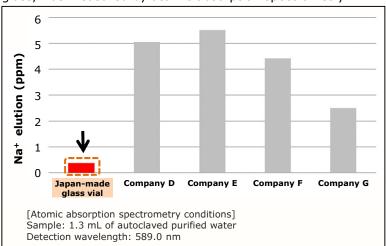

Japan-made glass vial was shown to be associated with low Na⁺ elution. As elution of alkaline components from vials causes sample decomposition, etc., it is critical to pay attention not only to sample adsorption to vials, but also elution from vials.

(3) Comparison of Impurities Elution from Septum Caps

Impurities elution from septum caps was tested to compare our Japan-made septum cap with septum cap products from other companies. Leachables in each vial were measured by GC/MS 10 times consecutively.

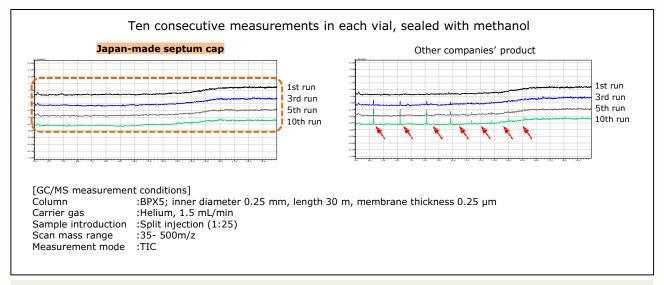

In consecutive measurements, leachables observed with other companies' products were hardly detected with Japan-made septum cap. Leachables from a cap or septum can have unexpected effects on test results, such as appearance of ghost peaks. Thus, it is necessary in analytical testing to pay attention to the quality not only of vials, but also of septum caps.

Product Information

Manufacturers of compatible products: Shimadzu, Hitachi, Waters, Agilent, etc.

Glass material :Borosilicate glass *low adsorption, low Na+ elution

Septum material :Fluororesin/silicone *low impurities elution

Specification of screw-top :9-425

List of JQ Vial Kit Products

All JQ Vial Kit products contain a Japan-made 1.5 mL screw-top vial.

Code No.	Vial Material	Write-on Spot	Preslit	Cap Color	Units per Package
636-52611	Clear Glass	-	-	Green	100
636-52591	Clear Glass	-	-	Blue	100
630-52631	Clear Glass	-	-	Black	100
634-52651	Clear Glass	-	-	White	100
633-52621	Clear Glass	-	•	Green	100
639-52601	Clear Glass	-	•	Blue	100
637-52641	Clear Glass	-	•	Black	100
631-52661	Clear Glass	-	•	White	100

Vial Material	Write-on Spot	Pre-slit	Cap Color	Units per Package
Clear Glass	•	-	Green	100
Clear Glass	•	-	Blue	100
Clear Glass	•	-	Black	100
Clear Glass	•	-	O White	100
Clear Glass	•	•	Green	100
Clear Glass	•	•	Blue	100
Clear Glass	•	•	Black	100
Clear Glass	•	•	O White	100
Amber Glass	-	-	Green	100
Amber Glass	-	-	Blue	100
Amber Glass	-	-	Black	100
Amber Glass	-	-	O White	100
Amber Glass	-	•	Green	100
Amber Glass	-	•	Blue	100
Amber Glass	-	•	Black	100
Amber Glass	-	•	White	100
Amber Glass	•	-	Green	100
Amber Glass	•	-	Blue	100
Amber Glass	•	-	Black	100
Amber Glass	•	-	White	100
Amber Glass	•	•	Green	100
Amber Glass	•	•	Blue	100
Amber Glass	•	•	Black	100
Amber Glass	•	•	White	100
	Clear Glass Amber Glass	Clear Glass Amber Glass	Clear Glass Clear	Clear Glass - Green Clear Glass - Blue Clear Glass - Black Clear Glass - White Clear Glass - Green Clear Glass - Blue Clear Glass - Black Clear Glass - White Amber Glass - Green Amber Glass - Blue Amber Glass - White Amber Glass - Blue Amber Glass - Black Amber Glass - Blue Amber Glass - Blu

Listed products are intended for laboratory research use only, and not to be used for drug, food or human use. / Please visit FUJIFILM Wako Laboratory Chemicals site: https://labchem-wako.fujifilm.com/ / This leaflet may contain products that cannot be exported to your country due to regulations. / Bulk quote requests for some products are welcomed. Please contact us.

FUJIFILM Wako Laboratory Chemicals site https://labchem-wako.fujifilm.com

FUJIFILM Wako Pure Chemical Corporation 1-2, Doshomachi 3-Chome, Chuo-ku, Osaka 540-8605, Japan Tel: +81 6 6203 3741 Fax: +81 6 6203 1999 ffwk-cservise@fujifilm.com

FUJIFILM Wako Chemicals U.S.A. Corporation 1600 Bellwood Road, Richmond, VA 23237, U.S.A. Toll-Free (U.S. only): +1 877 714 1920 Tel: +1 804 271 7677 Fax: +1 804 271 7791 wkuslabchem@fujifilm.com

FUJIFILM Wako Chemicals (Hong Kong) Limited FUJIFILM Wako (Guangzhou) Trading Corporation Room 1111, 11/F, International Trade Centre, 11-19 Sha Tsui Road, Tsuen Wan, N.T., Hong Kong Tel: +852-2799-9019 Fax: +852-2799-9808 wkhk.info@fujifilm.com

FUJIFILM Wako Chemicals Europe GmbH

Fuggerstr 12, 41468 Neuss, Germany Tel: +49 2131 311 0 Fax: +49 2131 311 100 labchem_wkeu@fujifilm.com

wkgz.info@fujifilm.com

Room 3003, 30/F., Dong Shan Plaza 69, Xian Lie Zhong Road, Guangzhou, 510095, China Tel: +86-20-8732-6381(Guangzhou) Tel: +86-21-6288-4751(Shanghai) Tel: +86-10-6413-6388(Beijing)