




## iMatrix-51

Substrate for cell culture iMatrix-511 is recombinant Laminin511-E8 fragments.

## Instructions

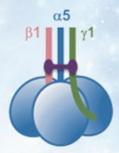

Laminin-511 E8 fragments

- 1) Dilute the solution with PBS (-). Coat dishes with 0.5 µg/cm<sup>2</sup>.
- 2) Allow the protein to bind for 3 h at room temperature.
- 3) Remove excess fluid from the coated surface.
- 4) Immediately plate the cells at desired density.

iMatrix-511 is a suitable substrate for ES/iPS cells. iMatrix-511 allows for the survival and xeno-free long-term single-cell passaging of ES/iPS cells.



## REFERENCES

Ido H et al. The requirement of the glutamic acid residue at the third position from the carboxyl termini of the laminin gamma chains in integrin binding by laminins. J. Biol. Chem. 282 (15): 11144-54, 2007

Taniguchi Y et al. The C-terminal region of laminin beta chains modulates the integrin binding affinities of laminins. J. Biol. Chem. 284 (12): 7820-31, 2009

Miyazaki T et al. Laminin E8 fragments support efficient adhesion and expansion of dissociated human pluripotent stem cells. Nat. Commun. 3: 1236. 2012

Nakagawa M et al. A novel efficient feeder-free culture system for the derivation of human induced pluripotent stem cells. Sci. Rep. 4: 3594, 2014

Doi D et al. Isolation of Human Induced Pluripotent Stem Cell-Derived Dopaminergic Progenitors by Cell Sorting for Successful Transplantation. Stem Cell Reports. 2 (3): 337-50, 2014

Takashima Y et al. Resetting Transcription Factor Control Circuitry toward Ground-State Pluripotency in Human. Cell. 158 (6): 1254-69, 2014

Fukuta M et al. Derivation of mesenchymal stromal cells from pluripotent stem cells through a neural crest lineage using small molecule compounds with defined media. PLos One. 9 (12): e112291, 2014

Burridge PW et al. Chemically defined generation of human cardiomyocytes. Nat. Methods. 11: 855-60, 2014

Okumura N et al. Laminin-511 and -521 Enable Efficient In Vitro Expansion of Human Corneal Endothelial Cells. Invest Ophthalmol Vis Sci. 56 (5), 2933-42, 2015

Sasaki K et al. Robust In Vitro Induction of Human Germ Cell Fate from Pluripotent Stem Cells. Cell Stem Cell. 17 (2), 178-94, 2015

Hayashi R et al. Co-ordinated ocular development from human iPS cells and recovery of corneal function. *Nature* 531, 367-80, 2016

Matsuno K et al., Redefining definitive endoderm subtypes by robust induction of human induced pluripotent stemcells Differentiation 2016.04.002

Nishimura K et al., Estradiol Facilitates Functional Integration of iPSC-Derived Dopaminergic Neurons into Striatal Neuronal Circuits via Activation of Integrin a5b1, Stem Cell Reports 6 (4) 511–524, 2016

Takayama K et al., Laminin 411 and 511 promote the cholangiocyte differentiation of human induced pluripotent stem cells, Biochemical and Biophysical Research Communications 474 (1): 91-96, 2016

	Product No.	Contents
iMatrix-511	892 011	350 $\mu$ g (500 $\mu$ g/mL) : 175 $\mu$ g×2 pcs.
	892 012	1,050 $\mu$ g (500 $\mu$ g/mL): 175 $\mu$ g×6 pcs.

You can coat more than 35 6-well plates with 1 mg iMatrix-511 for normal use.

Manufactured by

## Nippi, incorporated

1-1-1 Senjumidori-cho, Adachi-ku, Tokyo 120-8601, JAPAN PHONE: +81-3-3888-5184. FAX: +81-3-3888-5136 E-mail: protein-info@nippi-inc.co.jp

Distributor MATRIXOME, Inc.

3-2 Yamadaoka, Suita-shi, Osaka 565-0871, JAPAN Institute for Protein Research, Osaka University PHONE: +81-6-6877-0222. E-mail: info@matrixome.co.jp