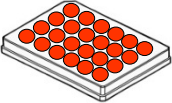
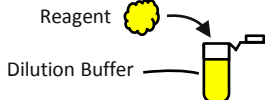
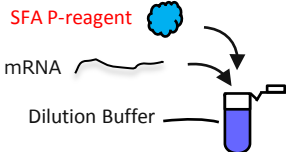
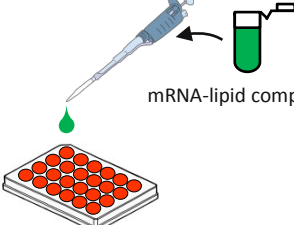
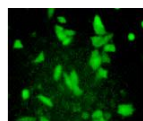
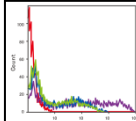


ScreenFect™A *plus* Transfection Protocol

The optimal condition for successful transfection varies. Start any new transfection by testing the recommended two concentrations of ScreenFectTM *plus* Reagent to determine an optimum amount.

2-Step method (Forward transfection method)

Timeline		Steps	Procedure Details																																																																
1	Day 0	 Pre-Cultured cells	Seed cells to be 70-90% confluent at transfection	<table><tr><th>Component</th><th colspan="2">96-well</th><th colspan="2">24-well</th><th colspan="2">12-well</th><th colspan="2">6-well</th></tr><tr><td>Adherent cells or suspension cells</td><td colspan="2">1.0-4.0 × 10⁴</td><td colspan="2">0.5-2.0 × 10⁵</td><td colspan="2">1.0-4.0 × 10⁵</td><td colspan="2">0.25-1.0 × 10⁶</td></tr></table> Seed cells to be 70-90% confluent at transfection. The medium replacement may improve a transfection efficiency in a 2-Step method.										Component	96-well		24-well		12-well		6-well		Adherent cells or suspension cells	1.0-4.0 × 10 ⁴		0.5-2.0 × 10 ⁵		1.0-4.0 × 10 ⁵		0.25-1.0 × 10 ⁶																																					
		Component	96-well		24-well		12-well		6-well																																																										
Adherent cells or suspension cells	1.0-4.0 × 10 ⁴		0.5-2.0 × 10 ⁵		1.0-4.0 × 10 ⁵		0.25-1.0 × 10 ⁶																																																												
2	Day 1	 Reagent Dilution Buffer	Dilute ScreenFect™A plus Reagent※ ¹ in Dilution Buffer, and then mix well ※ ¹ Vortex the reagent before use	<table><tr><th>Dilution Buffer for ScreenFect™A plus</th><th colspan="2">5 μL</th><th colspan="2">25 μL</th><th colspan="2">50 μL</th><th colspan="2">125 μL</th></tr><tr><td>mRNA : Transfection Reagent ratio</td><td>1 : 3</td><td>1 : 4</td><td>1 : 3</td><td>1 : 4</td><td>1 : 3</td><td>1 : 4</td><td>1 : 3</td><td>1 : 4</td></tr><tr><td>ScreenFect™A plus Transfection Reagent</td><td>0.15 μL</td><td>0.2 μL</td><td>0.75 μL</td><td>1.0 μL</td><td>1.5 μL</td><td>2.0 μL</td><td>3.75 μL</td><td>5.0 μL</td></tr></table>										Dilution Buffer for ScreenFect™A plus	5 μL		25 μL		50 μL		125 μL		mRNA : Transfection Reagent ratio	1 : 3	1 : 4	1 : 3	1 : 4	1 : 3	1 : 4	1 : 3	1 : 4	ScreenFect™A plus Transfection Reagent	0.15 μL	0.2 μL	0.75 μL	1.0 μL	1.5 μL	2.0 μL	3.75 μL	5.0 μL																											
		Dilution Buffer for ScreenFect™A plus	5 μL		25 μL		50 μL		125 μL																																																										
mRNA : Transfection Reagent ratio	1 : 3	1 : 4	1 : 3	1 : 4	1 : 3	1 : 4	1 : 3	1 : 4																																																											
ScreenFect™A plus Transfection Reagent	0.15 μL	0.2 μL	0.75 μL	1.0 μL	1.5 μL	2.0 μL	3.75 μL	5.0 μL																																																											
3	Day 1	 SFA P-reagent mRNA Dilution Buffer	Dilute mRNA and SFA P-reagent in Dilution Buffer, and then mix well	<table><tr><th>Dilution Buffer for ScreenFect™A plus</th><th colspan="2">5 μL</th><th colspan="2">25 μL</th><th colspan="2">50 μL</th><th colspan="2">125 μL</th></tr><tr><td>mRNA (0.1-2.5 μg / μL)</td><td colspan="2">50 ng</td><td colspan="2">250 ng</td><td colspan="2">500 ng</td><td colspan="2">1250 ng</td></tr><tr><td>SFA P-reagent (2μL / μg mRNA)</td><td colspan="2">0.1 μL</td><td colspan="2">0.5 μL</td><td colspan="2">1.0 μL</td><td colspan="2">2.5 μL</td></tr></table>										Dilution Buffer for ScreenFect™A plus	5 μL		25 μL		50 μL		125 μL		mRNA (0.1-2.5 μg / μL)	50 ng		250 ng		500 ng		1250 ng		SFA P-reagent (2μL / μg mRNA)	0.1 μL		0.5 μL		1.0 μL		2.5 μL																												
		Dilution Buffer for ScreenFect™A plus	5 μL		25 μL		50 μL		125 μL																																																										
mRNA (0.1-2.5 μg / μL)	50 ng		250 ng		500 ng		1250 ng																																																												
SFA P-reagent (2μL / μg mRNA)	0.1 μL		0.5 μL		1.0 μL		2.5 μL																																																												
4	Day 2	 mRNA-lipid complex	Add diluted mRNA (+SFA P-reagent) to diluted ScreenFect™A plus Reagent, and then incubate for 5 minutes ~ at room temperature※ ² ※ ² Incubation is available until the step 4 has been completed	<table><tr><td>Diluted mRNA (+SFA P-reagent)</td><td colspan="2">5 μL</td><td colspan="2">25 μL</td><td colspan="2">50 μL</td><td colspan="2">125 μL</td></tr><tr><td>Diluted ScreenFect™A plus Transfection Reagent</td><td colspan="2">5 μL</td><td colspan="2">25 μL</td><td colspan="2">50 μL</td><td colspan="2">125 μL</td></tr></table>										Diluted mRNA (+SFA P-reagent)	5 μL		25 μL		50 μL		125 μL		Diluted ScreenFect™A plus Transfection Reagent	5 μL		25 μL		50 μL		125 μL																																					
		Diluted mRNA (+SFA P-reagent)	5 μL		25 μL		50 μL		125 μL																																																										
Diluted ScreenFect™A plus Transfection Reagent	5 μL		25 μL		50 μL		125 μL																																																												
5	Day 2~	 mRNA-lipid complex	Add mRNA-lipid complex from step 3 to well of cell culture plate from step 1	<table><tr><th>Final composition [per well]</th><th colspan="2">96-well</th><th colspan="2">24-well</th><th colspan="2">12-well</th><th colspan="2">6-well</th></tr><tr><td>mRNA-lipid complex</td><td colspan="2">10 μL</td><td colspan="2">50 μL</td><td colspan="2">100 μL</td><td colspan="2">250 μL</td></tr><tr><td>mRNA amount</td><td colspan="2">50 ng</td><td colspan="2">250 ng</td><td colspan="2">500 ng</td><td colspan="2">1250 ng</td></tr><tr><td>SFA P-reagent used</td><td colspan="2">0.1 μL</td><td colspan="2">0.5 μL</td><td colspan="2">1.0 μL</td><td colspan="2">2.5 μL</td></tr><tr><td>ScreenFect™A plus Transfection Reagent used</td><td colspan="2">0.15 or 0.2 μL</td><td colspan="2">0.75 or 1.0 μL</td><td colspan="2">1.5 or 2.0 μL</td><td colspan="2">3.75 or 5.0 μL</td></tr><tr><td>Medium volume</td><td colspan="2">100 μL</td><td colspan="2">500 μL</td><td colspan="2">1000 μL</td><td colspan="2">2000 μL</td></tr></table>										Final composition [per well]	96-well		24-well		12-well		6-well		mRNA-lipid complex	10 μL		50 μL		100 μL		250 μL		mRNA amount	50 ng		250 ng		500 ng		1250 ng		SFA P-reagent used	0.1 μL		0.5 μL		1.0 μL		2.5 μL		ScreenFect™A plus Transfection Reagent used	0.15 or 0.2 μL		0.75 or 1.0 μL		1.5 or 2.0 μL		3.75 or 5.0 μL		Medium volume	100 μL		500 μL		1000 μL		2000 μL	
		Final composition [per well]	96-well		24-well		12-well		6-well																																																										
mRNA-lipid complex	10 μL		50 μL		100 μL		250 μL																																																												
mRNA amount	50 ng		250 ng		500 ng		1250 ng																																																												
SFA P-reagent used	0.1 μL		0.5 μL		1.0 μL		2.5 μL																																																												
ScreenFect™A plus Transfection Reagent used	0.15 or 0.2 μL		0.75 or 1.0 μL		1.5 or 2.0 μL		3.75 or 5.0 μL																																																												
Medium volume	100 μL		500 μL		1000 μL		2000 μL																																																												
		 	Visualize/analyze transfected cells	Incubate cells for 1 day ~ at 37°C. Then, analyze transfected cells.																																																															

For support, please visit the <http://screenfect.jp>